

Installing the Venable Matlab files

Move the Venload.m, Vendlmread.m and Vensave.m files to a directory in the Matlab path or create a directory and place it in the Matlab path. These files are installed in “C:\Program Files\Venable Instruments\Venable Software\Matlab m-files” for 32 bit Windows and they are installed in “C:\Program Files (x86)\Venable Instruments\Venable Software\Matlab m-files” for 64 bit Windows. In Windows 7, normal users cannot enter into Program Files. You can open a file explorer window to this directory from the Venable Software’s menu item “Help>>Matlab m-files”.

1. Add the directory to the Matlab path. You can do this by opening Matlab. For Matlab 5 and 6, select Set Path from the File menu. The Path Browser will open and you can add any directory to the Matlab path.
2. Make the “C:\Program Files\Venable Instruments\Venable Software\Matlab m-files” directory the current directory. For Matlab 5, use the CD command to change the current directory i.e. `cd 'C:\Program Files\Venable Instruments\Venable Software\Matlab m-files'`. For Matlab 6, select Current Directory from the View menu. The Current Directory Browser will open and you can make any directory the current directory.